


## ターゲットバージョン指定/解除 設定手順 (JP)

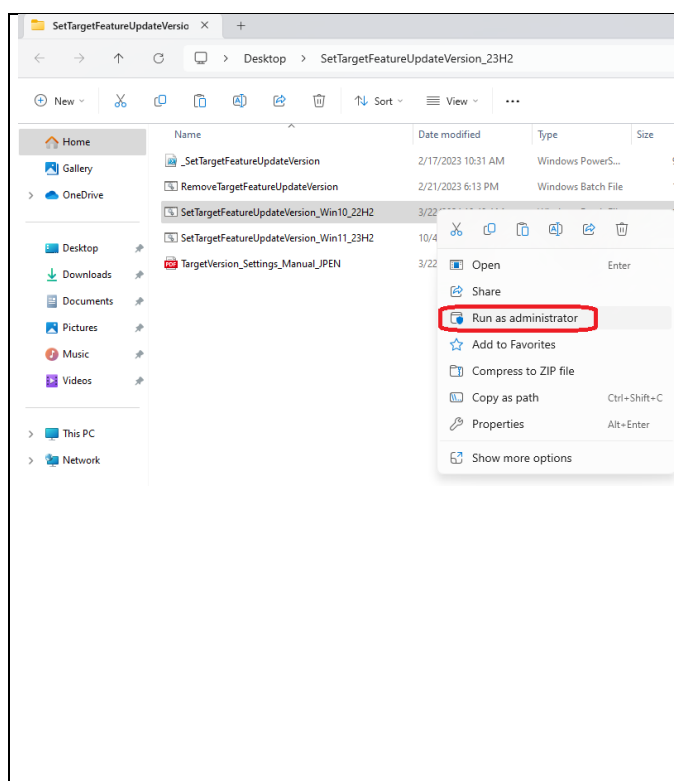

### 手順（１）

バッチファイルを右クリックし、表示されるコンテキストメニューからRun as Administratorを選択する。使用するバッチファイル名は以下の通りです。

ターゲットバージョン指定(Windows 10)  
用：


SetTargetFeatureUpdateVersion\_Win10\_22H2.bat

ターゲットバージョン指定(Windows 11)  
用：

SetTargetFeatureUpdateVersion\_Win11\_23H2.bat

ターゲットバージョン解除用：

RemoveTargetFeatureUpdateVersion.bat

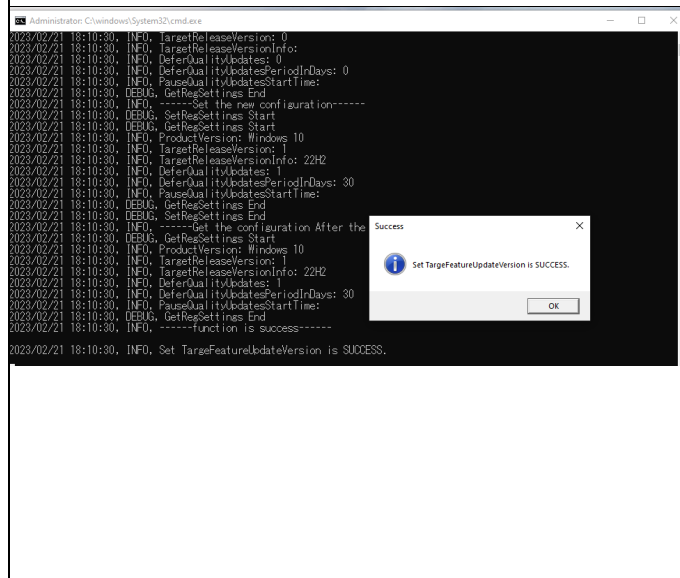



### 手順（２）

ダイアログ上で“SUCCESS”と表示されることを確認する。確認後、OKボタンでコマンドプロンプトを閉じて終了させる。

### （注１）

本バッチファイルと共通の設定をする [Local Group Policy Editor]内の[Select the target Feature Update version]設定がありますが、本バッチファイルで指定・解除を実行しても表示上の設定は変更されません。

## Set / Cancel Target version setting procedure (EN)

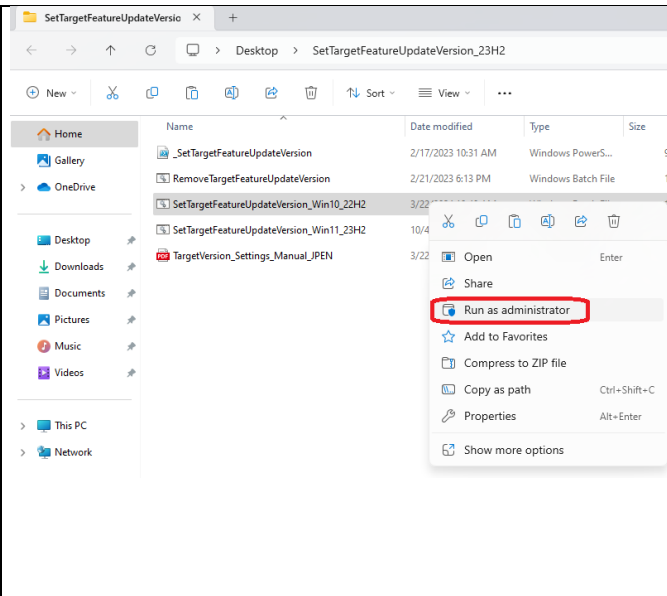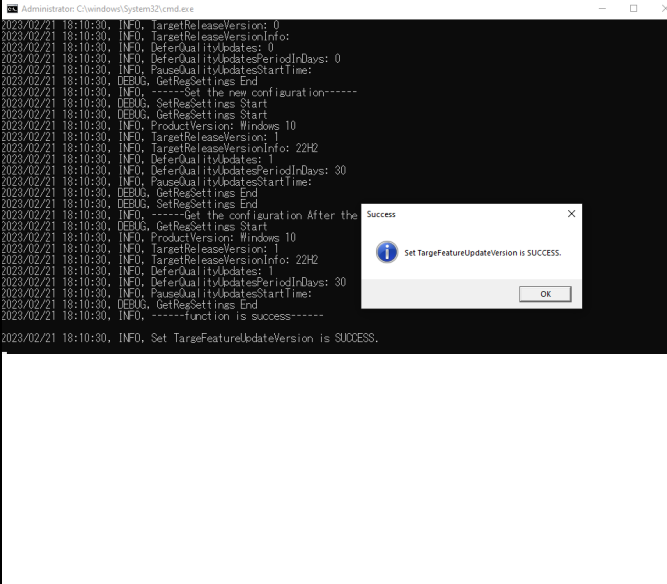	<p><b>Step 1</b></p> <p>Right-click the batch file and select “Run as Administrator” from the displayed context menu. The batch file name to be used is as follows.</p> <p>For target version setting (Windows 10): SetTargetFeatureUpdateVersion_Win10_22H2.bat</p> <p>For target version setting (Windows 11): SetTargetFeatureUpdateVersion_Win11_23H2.bat</p> <p>For canceling the target version setting: RemoveTargetFeatureUpdateVersion.bat</p>
	<p><b>Step 2</b></p> <p>Confirm that "SUCCESS" is displayed on the dialog. After confirmation, close the command prompt with the OK button and exit.</p> <p><b>Note:</b></p> <p>There is a [Select the target Feature Update version] setting in the [Local Group Policy Editor] that makes common settings with this batch file, but the settings on the display will not be changed even if you set / cancel in this batch file.</p>