

USB ライセンスキードライバインストール手順

注記 : Windows に管理者権限でログオンしてください。

注記 : 作業開始前に USB ライセンスキーが接続されていないことを確認して下さい。


注記 : 作業開始前に実行中のアプリケーションを全て終了してください。

1. ダウンロードしたファイル HASP_driver_v8_11.zip(または HASP_driver_v8_11)をインストールする環境のデスクトップにコピーします。
2. 手順 1 でコピーした圧縮ファイルを右クリックします。表示されたメニューから「すべて展開」を選択します。
圧縮ファイルは、展開すると以下のフォルダ構成になります。


-[展開先フォルダ]

- data
- Install.vbs (または Install)

3. 展開先フォルダの Install.vbs(または Install)をダブルクリックします。
4. 確認画面が表示されたら、「はい」を選択します。

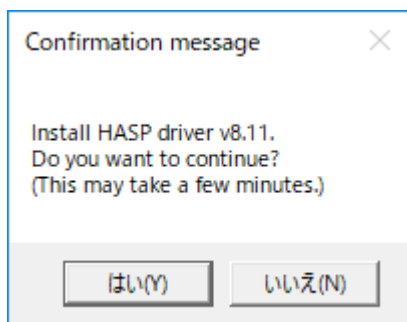


確認メッセージ

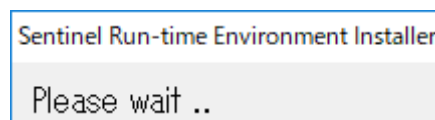

インストール画面

5. インストール完了画面が表示されたら、[OK]を選択します。USB ライセンスキードライバのインストールは完了です。

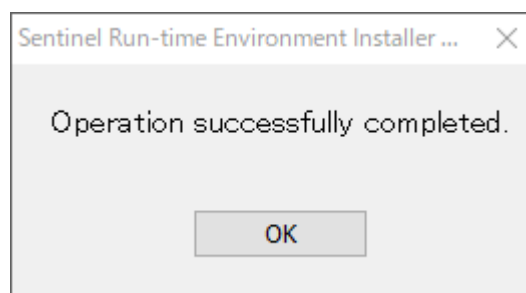

インストール完了画面

6. 手順 1 でデスクトップ上にコピーした修正パッチ圧縮ファイル及び、手順 2 で展開したフォルダ/ファイルを全て削除します。

以上


USB license key driver Installation Manual

NOTE : Log on to Windows in administrator privileges.


NOTE : Make sure that the USB license Key is not connected before starting work.

NOTE : Close all running applications before starting work.

1. Copy the compressed patch file "HASP_driver_v8_11.zip (or HASP_driver_v8_11)" to the Desktop of the environment to install USB license key driver.
2. Right-click the compressed patch file in the Step.1. Select [Extract All] from the menu displayed. The structure of the file after extraction is as follows.
 - [Extracted folder]
 - data
 - Install.vbs (or Install)
3. Double-click Install.vbs (or Install) of the extracted folder.
4. When the confirmation message appears, Select "Yes".

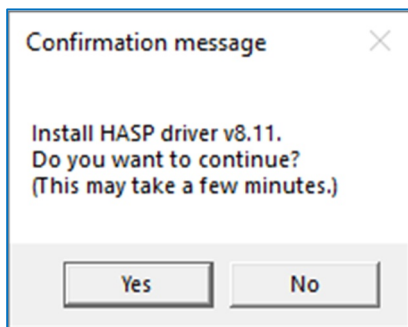


Confirmation message

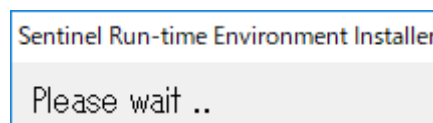

Install screen

5. When the installation completion screen appears, select "OK". The installation of USB license key driver is completed.

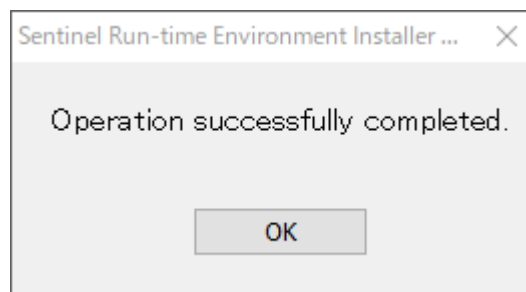

Installation completion screen

6. Delete all folders and files in the Step.1 and Step.2 on the Desktop.

End